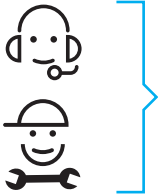
	EN Introduction DE Einleitung ES Introducción	FR Introduction ZH 简介 JA はじめに..... 4
	EN Safety Information DE Sicherheitshinweise ES Información de seguridad	FR Informations de sécurité ZH 安全信息 JA 安全情報..... 7
	EN Scope of delivery DE Lieferumfang ES Contenido de la entrega	FR Contenu de la livraison ZH 交货清单 JA 標準付属品..... 9
	EN Commissioning DE Inbetriebnahme ES Puesta en marcha	FR Mise en service ZH 调试 JA 試運転..... 10
	EN Pipette Check DE Pipettentest ES Comprobación de pipetas	FR Vérification des pipettes ZH 移液器检查 JA ピペット点検..... 13
	EN Maintenance DE Wartung ES Mantenimiento	FR Maintenance ZH 维护 JA メンテナンス..... 16
	EN Function test DE Funktionstest ES Prueba de funcionamiento	FR Test de fonctionnalité ZH 功能测试 JA 機能テスト..... 17
	EN Accessories and Spare Parts DE Zubehör und Ersatzteile ES Accesorios y piezas de	FR Accessoires et pièces de ZH 配件及备件 JA アクセサリおよびスペアパーツ..... 20
	EN Disposal DE Entsorgung ES Eliminación de residuos	FR Mise au rebut ZH 处置 JA 廃棄..... 22
	EN Technical data DE Technische Daten ES Características técnicas	FR Caractéristiques techniques ZH 技术资料 JA 技術データ..... 23
	EN Dimensions DE Abmessungen ES Dimensiones	FR Dimensions ZH 外形尺寸 JA 寸法..... 24
	EN Compliance information DE Informationen zur Konformität ES Información sobre conformidad	FR Information concernant la conformité ZH 合规性信息 JA コンプライアンス情報..... 25
	EN Warranty DE Gewährleistung ES Garantía	FR Garantie ZH 保修 JA 保証..... 26

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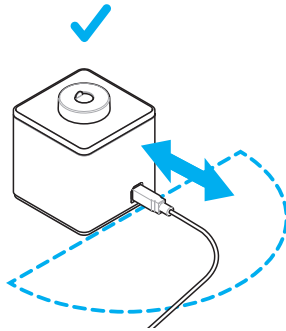
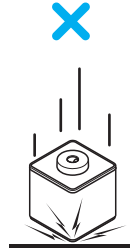
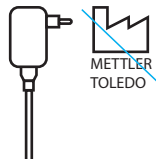
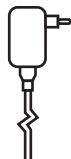
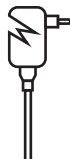
?

EN Intended use

SmartCheck™ is a pipette check instrument. It may be used for single channel or multichannel pipettes. The intended use is to check if a pipette's delivered volume meets the set instrument tolerances. SmartCheck™ shall be used for pipette checks only and cannot be used for pipette calibration. It is only to be used with distilled water. Any other type of use and operation beyond the limits of use stated by Mettler-Toledo GmbH without consent from Mettler-Toledo GmbH is considered as not intended.

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National approval documents, e.g., the FCC Supplier Declaration of Conformity, are available online and/or included in the packaging.

► <http://www.mt.com/ComplianceSearch>

Contact METTLER TOLEDO for questions about the country-specific compliance of your instrument.

United States of America

This equipment has been tested and found to comply with the limits for a Class A digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference when the equipment is operated in a commercial environment. This equipment generates, uses, and can radiate radio frequency energy and, if not installed and used in accordance with the instruction manual, may cause harmful interference to radio communications. Operation of this equipment in a residential area is likely to cause harmful interference in which case the user will be required to correct the interference at his own expense.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

1. This device may not cause harmful interference.
2. This device must accept any interference received, including interference that may cause undesired operation.

Changes or modifications not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

The FCC Supplier Declaration of Conformity is available online.

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FR Garantie
ZH 保修
JA 保証

EN Limited warranty and limitation of liability

Seller warrants the merchandise to conform to factory specifications. As buyer's exclusive remedy in the event of a breach of warranty, seller shall repair or replace, at its option, any nonconforming merchandise or parts there of for a period on one (1) year after delivery. For more info, visit www.mt.com/rainin

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GPP®

Good Pipetting Practice™

GPP® is a global pipette standard, ensuring consistent accuracy of pipetting processes, applicable to any pipette.

It helps to:

- Choose the appropriate pipette and tips
- Gain the range of pipetting skills necessary to produce reliable data
- Recognize the risk associated with out-of-calibration pipettes and the role of routine pipette checks vs. professional pipette service

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© Mettler-Toledo GmbH 03/2020